

FOUNDRY-MASTER UV

Metal Analysis – fast, easy and precise

The **FOUNDRY-MASTER** is a reliable, precise laboratory spectrometer for the qualitative and quantitative element analysis of metallic samples. The instrument is designed for stationary use as a bench-top-unit. The instrument is well established with hundreds of customers worldwide opting for its high performance qualities and practical features.

The instrument is based Optical Emission Spectroscopy (OES), the analysing method favoured by most metal producing and processing companies. The digital source (spark generator) is controlled via the external Windows™ PC and offers ideal excitation parameters for the most diverse alloys.

The high-resolution Multi-CCD-Optic utilises a traditional, robust, vacuum technology chamber, rather than economical and eccentric ways of removing harmful atmosphere from the optic. The optic range covers the complete wavelength range of 160 nm to 800 nm.

The rugged and approved vacuum technology is essentially maintenance-free and has significantly lower operating costs than alternative inert gas purged systems. The sample stand, open on three sides, offers the ability to measure almost any size or shape of sample, particularly large sample geometries. The unique design needs less cleaning than comparable systems, giving much longer operational time.

The cleaning process is fast and instrument recovery efficient due to pre-aligned parts and quick connectors. The **FOUNDRY-MASTER** is equipped with the unique WAS Jet-Stream Technology within the sample stand. The electrode is encapsulated in a flow of argon, providing two key benefits compared to previously available systems.

Argon consumption is reduced to a minimum, saving a considerable amount of operating expenses. Additionally, the sample does not need to completely seal the spark stand opening. Tubes, bars and even wires can be directly analysed, using only one universal adapter. This is a major operational advantage and strongly reduces the sample preparation time.

The **FOUNDRY-MASTER** can be upgraded with new calibration modules at any time and without expensive and time-consuming modifications. Low argon consumption thanks to optimisation of argon flush and use of vacuum technology. Analysis of all common metals and their alloys.

Three side open spark stand for measuring large sample geometries. Jet-Stream Technology for the precise analysis of curved and irregular surfaces, including wires.

- Analysis of all metals and alloys
- Jet-Stream Technology
- Low argon consumption
- Precise analysis including material identification
- Easy-access spark stand



The Business of Science®

Worldwide Analytical Systems AG



an Oxford Instruments company

OXFORD
INSTRUMENTS

- Analysis of all metals and alloys
- Jet-Stream Technology
- Low argon consumption
- Precise analysis including material identification
- Easy-access spark stand

Oxford Instruments Industrial Analysis

UK

Halifax Road, High Wycombe
Bucks, HP12 3SE England
Tel: +44 (0) 1494 442255
Fax: +44 (0) 1494 461033
Email: analytical@oxinst.com

China

Beijing
Tel: (8610) 6518 8160/1/2
Fax: (8610) 6518 8155
Email: info@oichina.cn

Finland

Espoo
Tel: +358 9 329 411
Fax: +358 9 3294 1300
Email: FI-Espoo_Info@oxinst.com

France

Saclay, Cedex
Tel: +33 (0) 1 69 85 25 24
Fax: +33 (0) 1 69 41 86 80
Email: analytical-info@oxford-instruments.fr

Germany

WAS Worldwide Analytical Systems AG
Wellesweg 31 D-47589 Uedem
Tel: +49 (0) 2825 9383 0
Fax: +49 (0) 2825 9383 100
Email: info@was-ag.com
www.was-ag.com

Japan

Tokyo
Tel: +81 (0) 3 5245 3591
Fax: +81 (0) 3 5245 4466/4477
Email: oikkma@oxinst.co.jp

Latin America

Clearwater FL
Tel: +1 727 538 7702
Fax +1 727 538 4205
Email: oxford@gate.net

Singapore

Tel: +65 6337 6848
Fax: +65 6337 6286
Email: xrf.sales@oxford-instruments.com.sg

North America

Elk Grove Village IL
Tel: +1 847 439 4404
Fax: +1 847 439 4425
Email: sales@msys.oxinst.com

www.oxford-instruments.com



an Oxford Instruments company



Technical Data

Height	368 mm (14,5")
Width	889 mm (35,0")
Depth	635 mm (25,0")
Weight	70 kg (154,3 lbs)
Mains Power	230-110 VAC (50/60 Hz)
Operating Mode	600 W
Standby Mode	50 W

Solid State Source

Computer controlled parameters	
Frequency	100 – 400 Hz
Voltage	300 – 500 V
High Energy Pre Spark HEPS	

Readout System

External PC-Workstation incorporating up-to-date technology

Vacuum System

Direct-path optical spark-stand
Low-noise vacuum pump
Shut-off-valve for easy replacement of the optical window

Options

Wire-adaptor set
Sample preparation devices
Spare parts kit
Consumables kit

Optical System

Multi CCD optical-system with Paschen-runge Mount	
Resolution CCD	6 Pico-meter
Reciprocal dispersion	0,9 nm/mm (1st order)
Focal length	350 mm
Holographic grating	3000 g/mm
Wavelength	160 – 800 nm

Typical Applications

Analytical mode / identification
Majority of metals and their alloys
Fe- alloys Cast-iron alloys
Al: alloys, cast alloys...
Cu: bronze, brass, Cu – Ni,...
Ni: hastelloy ~ inconel ~ monel,...
Ti: ti pure . Ti.6 - 4 ~ Ti.8-Mn...
Mg-, Zn - alloys, solders and many more...

We put the spark in Spectrometry

Click onto www.oxford-instruments.com for more information

Oxford Instruments, at High Wycombe, UK, operates Quality Management Systems approved to the requirements of BS EN ISO 9001. This publication is the copyright of Oxford Instruments Analytical Limited and provides outline information only which (unless agreed by the company in writing) may not be used, applied or reproduced for any purpose or form part of any order or contract or be regarded as a representation relating to the products or services concerned. Oxford Instruments' policy is one of continued improvement. The company reserves the right to alter, without notice, the specification, design or conditions of supply of any product or service. Oxford Instruments acknowledges all trade marks and registrations.

© Oxford Instruments Analytical Ltd, 2008. All rights reserved.

As part of Oxford Instruments' environmental policy this brochure has been printed on FSC paper.



Certificate No FM29142

Part no: OIIA/029/A/0108